



"Standards for the Instrument Shop and the Metrology Lab"

For a detailed listing of our products, with full specification sheets in pdf format, please visit our website at www.calibrationsales.com Over 5000 products for the calibration lab.

62 Cove Road, Stonington, CT 06378

Tel: 860/572-0111 Fax: 860/572-7744 Email: dan@calibrationsales.com

Thunder Scientific Model 2500

Mobile "Two-Pressure" Humidity Generator

FEATURES

- Accuracy ± 0.5 %RH
- Traceable to NIST
- Self Contained and Mobile
- Automated Control of User Setpoints
- Shipped with ControLog™ Automation Software and HumiCalc® Humidity Conversion Software
- Computerized Internal Transducer Calibration
- Automatically Applies Enhancement Factors
- RS-232C Serial Interface



DESCRIPTION

The **Model 2500** Humidity Generator is a self contained system capable of producing atmospheres of known humidities using the fundamental, NIST proven, "two-pressure" principle. This system is capable of continuously supplying accurately known humidity values for instrument calibration, evaluation, and verification, as well as for environmental testing.

Simply apply power and the **2500** will power-up ready to generate. Humidity setpoint values are input by the operator from the front panel keypad and are limited only by the range of the **2500** humidity generator.

Relative humidities are calculated from the measurements of pressure and temperature with the following formula:

$$\text{Formula: \%RH} = f_s/f_c \times e_s/e_c \times P_c/P_s \times 100$$

To generate a known humidity, the computer controls the pressure ratio P_c/P_s utilizing the enhancement factor ratio f_s/f_c and the effective degree of saturation e_s/e_c . Humidity

			SetPnt	Actual	
% RH	@ Pc		20.05	20.05	CHNG SETP
*% RH	@ PcTc		20.00	20.00	CHNG UNIT
SATUR	PSI	€1.40	€1.40		EDIT /CAL
CHMBR	PSI			12.17	
SATUR	°C	23.00	23.00		
CHMBR	°C			23.05	
FLOW	l/m	20.00		13.88	RUN
09/09/99		11:45:18		■■■■■JF	

produced is solely dependent on the measurement of pressures and temperatures and does not rely on any other device (such as a dew point hygrometer, psychrometer, or humidity sensor) for the measurement of water vapor content. Precision humidity generation is determined by the accuracy of the pressure measurements and on the accuracy and uniformity of temperature throughout the generating system.

PRINCIPLE OF OPERATION

The **Model 2500** Humidity Generator operates using an on board multifunction CPU in conjunction with other peripheral cards to perform calculation and control functions. The embedded computer control system allows the **2500** to generate known humidity levels unattended, freeing the operating technician from the task of system monitoring and adjustment. A computer and/or printer may be connected via the bi-directional RS-232C interface ports allowing remote setpoint control and continuous acquisition of system data.

Humidity and temperature setpoint values are input by the operator from the front panel keypad while visual indications of system status are displayed in real time on the liquid crystal display.

Control / Display Screen


All Control and measurement parameters critical to the operation of the humidity generator are displayed on this screen. Each parameter in the leftmost column is identified with a brief title and corresponding units. The generator operates in a variety of user selectable pressure, temperature, and flow units. Some of these are °C, °F, psi, "Hg, Tor, mbar, kPa, l/m, l/h, cfm and cfh. Humidity is calculated and displayed in percent relative humidity (%RH). The asterisk in the leftmost column indicates the active humidity control parameter. The "SetPnt" column lists control setpoints and the "Actual" column lists all of the measured data and calculated parameters of the generator.

Temperature Control: The system utilizes a fluid jacketed test chamber for extremely stable temperature control. Temperature setpoint control is attained by controlling the temperature of the circulating fluid medium that jackets the test chamber and associated humidity generation components. Chamber and saturation temperatures are governed by this medium and are digitally controlled by the computer at any value between 0 °C and 70 °C using PID (proportional-integral-derivative) algorithms.

Pressure and Flow Control: Pressure control and mass flow control are accomplished through computer actuation of electromechanical valve assemblies. Pressure and flow are

measured continuously and controlled using PID algorithms similar to those employed in temperature control.

		Count	Deg C	
*Satur	Tmp	1180	23.41	LOW TEMP
*PrSat	Tmp	1165	23.12	MID TEMP
*Expan	Tmp	1173	23.20	HIGH TEMP
*Chamb	Tmp	1155	23.08	
Reter	Tmp	1257	.1364	
LOW		MID	HIGH	EXIT
0		35	70	QUIT

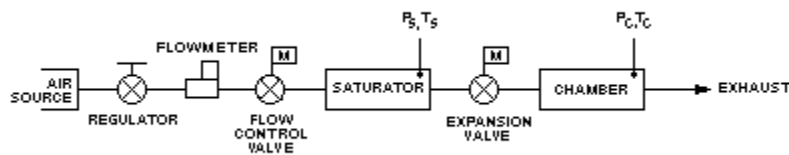

Elemental Schematic of the 2500 generator.

Calibration: The 2500 humidity generator employs an integral programmatic calibration scheme allowing the temperature and pressure transducers to be calibrated while they are electrically connected to the humidity generator.


Calibration / Display Screen

Coefficients for each transducer are calculated by the computer and stored in the system's nonvolatile memory until the next calibration is performed.



APPLICATIONS

The fluid jacketed test chamber can accommodate humidity sensors, hygrothermographs, chilled mirror hygrometers, and various material samples for environmental testing. Virtually any humidity and temperature point may be generated, for any length of time, within the operational limits of the generator. The output of the device under test may then be compared with the generator's printed data for analysis.

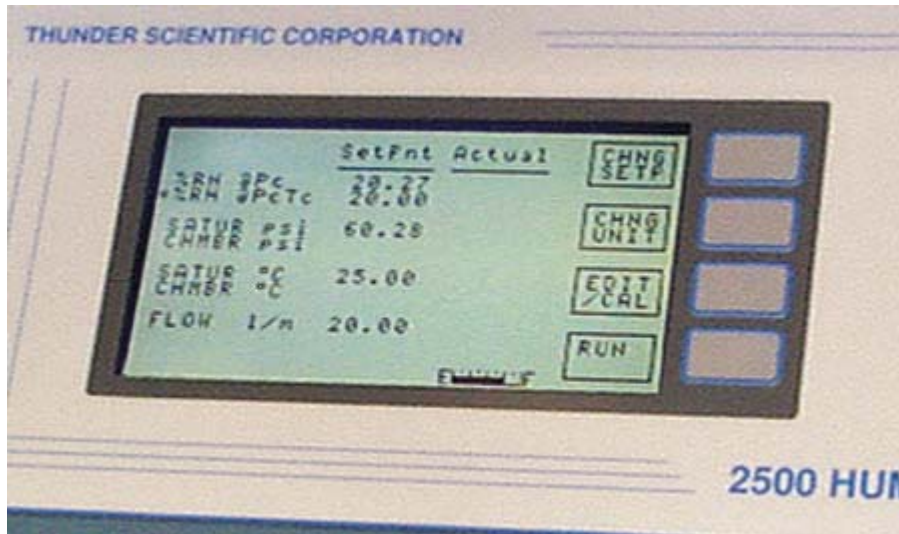

The liquid crystal display has many different user features, like the liquid level readout.

Humidity Sensors And Chart Recorders: Insert humidity probes through the two inch port in the side of the chamber or place hygrothermographs into the chamber and you can: determine humidity calibration accuracy and characterize humidity sensitivity by subjecting the sensing system to a variety of humidity levels; perform operational checks such as the sensing systems capability to correctly calculate and display other humidity parameters; determine the repeatability, stability, hysteresis, and drift characteristics of various humidity sensing systems.


Chilled Mirror Hygrometers: Install the actual chilled mirror head into the chamber or insert a sample tube through the test port and draw a sample through the chilled mirror head and you can: verify mirror temperature measurement accuracy (calibration) when the hygrometer is in thermal equilibrium with its environment; perform operational checks of the heatpump and optical components before and after mirror cleaning and balancing; determine whether the hygrometer is controlling the mirror deposit in the liquid phase or ice phase when operating at dew and frost points below 0 °C; determine if the hygrometer is correctly calculating other humidity parameters; determine the hygrometer's repeatability, stability, and drift characteristics.

Environmental Testing: The **2500** can serve as a test bed for evaluation and R&D of humidity sensors, humidity sensing systems, and humidity sensitive products, e.g., polymers, composites, film, magnetic medium, blood gas analysis, pharmaceuticals, soil hydrology, consumables, electronics, optics, etc. Depending on the temperature and humidity being generated, the **2500** may operate continuously from hours to months. With continuous generation of a nominal 50 %RH at 21 °C, the reservoir will last about two weeks between refills.

We've added a NEW model to our product line. The **Model ACS-517B** Oil-less Compressed Air System (For Use With The Model 2500).



The **Model ACS-517B** has a hose hold down on top for ease of storage.

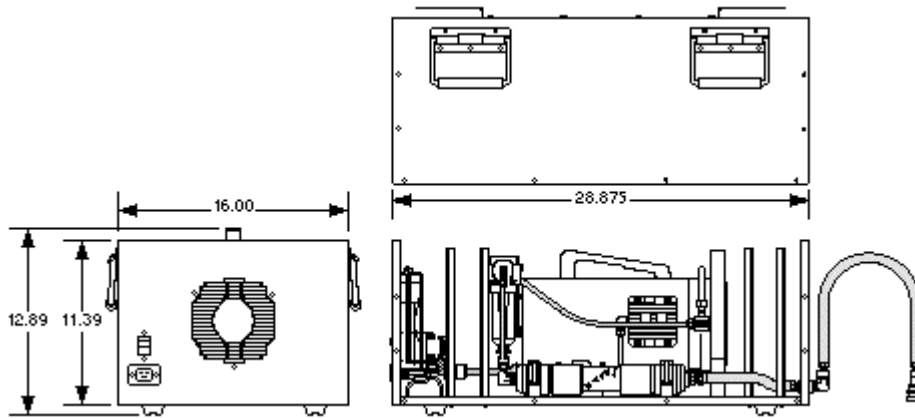

Physical Dimensions:

The **Model ACS-517B** is used with a Model 2500 and will fit on the bottom shelf of the 2500 mobile cart. This is a fully enclosed compressed air supply with an air dryer and sound muffling cabinet. This system is ideal for in Lab use because the sound level is less than 70 decibels.

Model 2500 Generator

SPECIFICATIONS

Relative Humidity Range: 10% to 98%
Relative Humidity Resolution: 0.02%
Relative Humidity Accuracy: $\pm 0.5\%$ @ $P_c T_c (k=2)$
Chamber Temperature Range: 0°C to 70°C
Chamber Temperature Range:(Optional) -10°C to +70°C
Chamber Temperature Resolution: 0.02°C
Chamber Temperature Uniformity:** $\pm 0.1^\circ\text{C}$
Chamber Temperature Accuracy: $\pm 0.06^\circ\text{C}$
Chamber Pressure Range: Ambient
Gas Flow Rate Range: 2 to 20 slpm
Gas Type: Air or Nitrogen
Gas Pressure Rating (MAWP): 175 psiG
Heating/Cooling Rate: 2.5 Minutes Per °C Average
Chamber Window: 6" x 6" (152 mm x 152 mm) Window in Door
Physical Dimensions: See Table A
Physical Dimensions of Cart: See Table B
Physical Dimensions 2500 With Cart: See Table C
Chamber Dimensions: See Table D
Access Port: See Table E

** When operated at temperatures within 10 °C of room ambient temperature.

See photos below...

TABLE A
Physical Dimensions

 <p align="center">MODEL</p>	H DIM	W DIM	D DIM
2500	19" (483 mm)	33" (838 mm)	20" (508 mm)
2500S	19" (483 mm)	36" (914 mm)	20" (508 mm)
2500ST	22" (559 mm)	36" (914 mm)	23" (584 mm)

Not including feet, handle, or other protrusions.



TABLE B
Physical Dimensions of Cart

MODEL	H DIM	W DIM	D DIM
--------------	--------------	--------------	--------------

2500	33.25" (845 mm)	40" (1.20 m)	23" (584 mm)
2500S	33.25" (845 mm)	43" (1.09 m)	23" (584 mm)
2500ST	33.25" (845 mm)	43" (1.09 m)	26" (660 mm)

TABLE C
Physical Dimensions 2500 with Cart

MODEL	H DIM	W DIM	D DIM
2500	53" (1.35 m)	40" (1.20 m)	23" (584 mm)
2500S	53" (1.35 m)	43" (1.09 m)	23" (584 mm)
2500ST	56" (1.42 m)	43" (1.09 m)	26" (660 mm)

TABLE D
Chamber Dimensions

MODEL	H DIM	W DIM	D DIM
2500	12" (305 mm)	12" (305 mm)	10" (254 mm)
2500S	12" (305 mm)	15" (381 mm)	10" (254 mm)
2500ST	15" (381 mm)	15" (381 mm)	12" (305 mm)

TABLE E
Access Port Dimensions

Option	# Ports	Port Diameter	Location
Standard	1	1.9" (48 mm)	Right Side
-TPA	2	1.9" (48 mm)	Right Side
	1	1/4" Swagelok	Right Side
-DPV	1	1.9" (48 mm)	Right Side
	1	1/4" Swagelok	Right Side
-MPD	6	1.1" (28 mm)	In Door

Other custom options are available.

UTILITIES

Electrical Power: 100/120 V~, 15 A, 50/60 Hz
(Optional) 200/240 V~, 8 A, 50/60 Hz
Air Compressor: 100/120 V~, 5 A, 50/60 Hz
(Optional) 200/240 V~, 2.5 A, 50/60 Hz
Air Supply (External): Clean, Oil Free Instrument Air
..... @ 175 psiG & 20 slpm

ENVIRONMENTAL

Operating Temperature: 15°C to 30°C
Storage Temperature: 0°C to 50°C
Humidity: 5 to 95% Non-condensing

For further information or a quote please contact:

HALL ASSOCIATES INC

62 COVE ROAD
STONINGTON, CT 06378
Attn: Dan Hall

TEL: 860/572-0111
FAX: 860/572-7744

[E-mail: dan@calibrationsales.com](mailto:dan@calibrationsales.com)